

---

## **CITRATE BUFFER (Ready-to-use) pH6.0**

### **(0.01 M Citrate Acid Buffer)**

Cat # NB-23-00172

Size: 1000ml liquid form.

#### **Recommended Protocol**

---

0.01 M Citrate Acid Buffer is ready-to-use. Refer to HIER Procedure for details.

#### **Precaution**

---

Wear gloves and take other necessary laboratory safety procedure.

#### **Storage**

---

Store at 2-8°C. Do not freeze.

#### **Procedure for FFPE Tissue Sections**

---

1. Deparaffinize in xylene and rehydrate tissue in graded alcohols.
2. Block endogenous peroxidase with hydrogen peroxide solution for 10 minutes.
3. Rinse with distilled water.
4. Rinse with PBS.
5. Heat ready-to-use citrate buffer in a beaker on a hot plate until the temperature of the buffer reaches 93°C-95°C. Place slides in heated buffer for 10-15 minutes.
6. Remove beaker with slides from the hot plate and allow it cool for 25-30 minutes.
7. Rinse slides with PBS 3 times, 2 minutes each time.
8. Apply avidin/biotin blocking or general blocking if required
9. Start immunostaining procedure.

*For reference only*

*For Research Use Only. Not for Diagnostic or Therapeutic Use.*