# cells in focus **bidi**®

## The ibidi Chambered Coverslips

Versatile µ-Slides for Advanced Microscopy


- ✓ Available with 1–18 wells for experimental flexibility
- Multiple surface options to suit your research needs
- ✓ Exceptional optical clarity for high-end imaging

"

l tested the μ-Slide 18 Well and the μ-Slide 8 Well<sup>high</sup>. I used them for confocal microscopy.

I can say that I am very pleased with these products! Super **easy to use** and **very convenient**.

> Rita Ribeiro University of Lisbon, Portugal

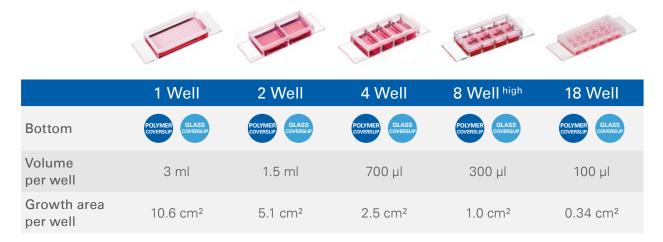

## The ibidi Chambered Coverslip Variety


# Distributed by: CliniSciences Group

## The ibidi Chambered Coverslips

Versatile µ-Slides for Advanced Microscopy

## The Principle of Imaging Chambers: The Coverslip Bottom

The outstanding characteristic of the ibidi  $\mu$ -Slides is their thin coverslip bottom, which has excellent features for high-end microscopy applications. In comparison, the bottom of standard cell culture plastics is about 1 mm thick—which is more than 5 times the thickness of the coverslip and, therefore, not ideal for imaging.

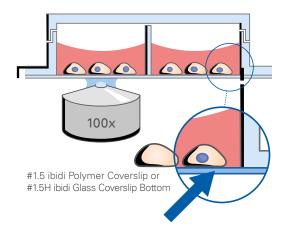



ibidi Polymer Coverslip


The ibidi Polymer Coverslip Bottom is suitable for various imaging techniques up to the highest resolution. With a standard #1.5 coverslip thickness of 180  $\mu$ m (+10/–5  $\mu$ m), it meets all optical requirements for microscopes. The ibidi Polymer Coverslip is compatible with a variety of immersion oils, which are specified at ibidi.com/oil.

# Surfaces and Coatings for the ibidi Polymer Coverslip




## ibiTreat (Tissue Culture-Treated)

Excellent adhesion of adherent cells, hydrophilic surface with no need for any additional coating.

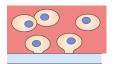

#### Hydrophobic, Uncoated Surface Weak adhesion of adherent cells, suitable for the application of specific coatings.

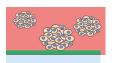

### Bioinert (ULA) Surface

No adhesion of adherent cells or any biomolecule, ideal for spheroid and organoid culture.

#### Coated Surface Culture of adhe Collagen I, IV,


Culture of adherent cells on a Collagen I, IV, or Poly-L-Lysine surface.

## ibidi Glass Coverslip




The ibidi Glass Coverslip Bottom was developed specifically for TIRF, super-resolution microscopy, and single molecule microscopy. However, it is also ideally suitable for standard imaging techniques. The D 263 M Schott borosilicate glass has a #1.5H thickness of 170  $\mu$ m (+/-5  $\mu$ m) and unrestricted immersion oil compatibility.

# Surfaces for the ibidi Glass Coverslip

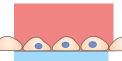

#### **Glass Surface**

Adhesion of adherent cells (coating might be required), ideal for special microscopy applications.

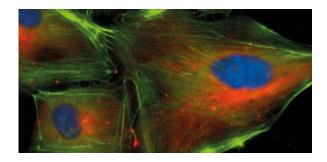

Immunofluorescence of human vascular endothelial cells (HUVECs) in a μ-Slide 18 Well ibiTreat. Red: alpha-Tubulin, green: F-actin, blue: nuclei.

# Distributed by: CliniSciences Group

# **CliniSciences** Group

#### Austria

Company: CliniSciences GmbH Address: Sternwartestrasse 76, A-1180 Wien - Austria Telephone: +43 720 115 580 Fax: +43 720 115 577 Email: oesterreich@clinisciences.com Web: https://www.clinisciences.com

#### Finland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064

Email: suomi@clinisciences.com Web: https://www.clinisciences.com

#### Iceland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: island@clinisciences.com Web: https://www.clinisciences.com

#### Netherlands

Company: CliniSciences B.V. Address: Gaetano Martinolaan 85, 6229 GS Maastricht - Netherlands Telephone: +31 85 2082 351 Fax: +31 85 2082 353 Email: nederland@clinisciences.com Web: https://www.clinisciences.com

#### Portugal

Company: Quimigen Unipessoal LDA Address: Rua Almada Negreiros, Lote 5, Loja 14, 2615-275 Alverca Do Ribatejo - Portugal Telephone: +351 30 8808 050 Fax: +351 30 8808 052 Email: info@quimigen.com Web: https://www.quimigen.pt

#### Switzerland

Company: CliniSciences AG Address: Fracht Ost Flughafen Kloten CH-8058 Zürich - Switzerland Telephone: +41 (044) 805 76 81 Fax: +41 (044) 805 76 75 Email: switzerland@clinisciences.com Web: https://www.clinisciences.com

#### Belgium

Company: CliniSciences S.R.L Address: Avenue Stalingrad 52, 1000 Brussels - Belgium Telephone: +32 2 31 50 800 Fax: +32 2 31 50 801 Email: belgium@clinisciences.com Web: https://www.clinisciences.com

### France

Company: CliniSciences S.A.S Address: 74 Rue des Suisses, 92000 Nanterre- France Telephone: +33 9 77 40 09 09 Fax: +33 9 77 40 10 11 Email: info@clinisciences.com Web: https://www.clinisciences.com

#### Ireland

Company: CliniSciences Limited Address: Ground Floor, 71 lower Baggot street Dublin D02 P593 - Ireland Telephone: +353 1 6971 146 Fax: +353 1 6971 147 Email: ireland@clinisciences.com Web: https://www.clinisciences.com

#### Norway


Company: CliniSciences AS Address: c/o MerVerdi Munkerudtunet 10 1164 Oslo - Norway Telephone: +47 21 988 882 Email: norge@clinisciences.com Web: https://www.clinisciences.com

Spain

Company: CliniSciences Lab Solutions Address: C/ Hermanos del Moral 13 (Bajo E), 28019, Madrid - Spain Telephone: +34 91 269 40 65 Fax: +34 91 269 40 74 Email: espana@clinisciences.com Web: https://www.clinisciences.com

#### υĸ

Company: CliniSciences Limited Address: 11 Progress Business center, Whittle Parkway, SL1 6DQ Slough- United Kingdom Telephone: +44 (0)1753 866 511 or +44 (0) 330 684 0982 Fax: +44 (0)1753 208 899 Email: uk@clinisciences.com IWeb: https://www.clinisciences.com

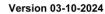

#### Denmark

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: danmark@clinisciences.com Web: https://www.clinisciences.com

#### Germany

Company: Biotrend Chemikalien GmbH Address: Wilhelm-Mauser-Str. 41-43, 50827 Köln - Germany Telephone: +49 221 9498 320 Fax: +49 221 9498 325 Email: info@biotrend.com Web: https://www.biotrend.com

#### Italy

Company: CliniSciences S.r.I Address: Via Maremmana inferiore 378 Roma 00012 Guidonia Montecelio - Italy Telephone: +39 06 94 80 56 71 Fax: +39 06 94 80 00 21 Email: italia@clinisciences.com Web: https://www.clinisciences.com

#### Poland

Company: CliniSciences sp.Z.o.o. Address: ul. Rotmistrza Witolda Pileckiego 67 lok. 200 - 02-781 Warszawa -Poland Telephone: +48 22 307 0535 Fax: +48 22 307 0532 Email: polska@clinisciences.com Web: https://www.clinisciences.com

#### Sweden

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: sverige@clinisciences.com Web: https://www.clinisciences.com

#### USA

Company: Biotrend Chemicals LLC Address: c/o Carr Riggs Ingram, 500 Grand Boulevard, Suite 210 Miramar Beach, FL 32550- USA Telephone: +1 850 650 7790 Fax: +1 850 650 4383 Email: info@biotrend-usa.com Web: https://www.biotrend-usa.com













